



Proficiency Test	Assigned value (mg/L)	Result (mg/L)	Z-score
BIPEA Food Allergens – EGG 77B - Pasta – 09 - 2577 - March 2024	NOT DETECTED	<0.4	-
BIPEA Food Allergens – EGG 77B - Pasta – 08 - 2677 - March 2024	22.2	24.2	0.29
BIPEA Food Allergens – EGG 77B - Pasta – 08 - 2777 - March 2024	54.1	63.6	0.74
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1077 – January 2024	NOT DETECTED	<0.4	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 16 - 1177 – January 2024	12.8	10.2	-0.51

BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 16 - 1277 - January 2024	17.0	14.3	-0.35
BIPEA Food Allergens - EGG 77B - Cake Mix - 09 - 0477 - November 2023	50.9	37.9	-0.71
BIPEA Food Allergens - EGG 77B - Cake Mix - 09 - 0577 - November 2023	17.0	16.2	-0.16
BIPEA Food Allergens - EGG 77B - Cake Mix - 09 - 0677 - November 2023	NOT DETECTED	<0.4	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 15 - 1077 - June 2023	NOT DETECTED	<0.4	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 15 - 1277 - June 2023	23.1	22.3	-0.09
BIPEA Food Allergens - EGG 77B - Pasta - 08 - 2577 - March 2023	40.5	50.8	0.68
BIPEA Food Allergens - EGG 77B - Pasta - 07 - 2677 - March 2023	NOT DETECTED	<0.4	-
BIPEA Food Allergens - EGG 77B - Pasta - 07 - 2777 - March 2023	19.5	27.6	0.94
BIPEA Allergens in Industrial Process Water 77D - 14 2022-2023 - 0381 Industrial Process Water - 14 - 1077 - January 2023	26.8	24.3	-0.23
BIPEA Allergens in Industrial Process Water 77D - 14 2022-2023 - 0381 Industrial Process Water - 14 - 1177 - January 2023	21.6	18.6	-0.35
BIPEA Allergens in Industrial Process Water 77D - 14 2022-2023 - 0381 Industrial Process Water - 14 - 1277 - January 2023	NOT DETECTED	<0.4	-

BIPEA Food Allergens – EGG 77B - Cake Mix – 08 – 0477 - October 2022	NOT DETECTED	<0.4	-
BIPEA Food Allergens – EGG 77B - Cake Mix – 08 – 0577 - October 2022	38.9	45.0	0.56
BIPEA Food Allergens – EGG 77B - Cake Mix – 08 – 0677 - October 2022	11.6	19.8	1.67
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 13 - 1077 - June 2022	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 13 - 1177 - June 2022	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water – 13 - 1277 - June 2022	NOT DETECTED	<0.4	-

<b>Proficiency Test</b>	<b>Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Egg Ei/Egg protein R6402</b>	<b>Range for <math> z  \leq 2</math> according to R-Biopharm Ridascreen Fast Egg Ei/Egg protein R6402</b>	<b>Result (mg/kg)</b>	<b>Recovery (%)</b>
FAPAS Allergens in Cooked Biscuit 27382 January – March 2024	61.5	30.7 – 92.2	47.6	77.4
FAPAS Allergens in Cake Mix - 27377A - November 2023 – January 2024	29.3	14.6 – 43.9	33.4	114.0
FAPAS Allergens in Cake Mix - 27377B - November 2023 – January 2024	36.1	18.0 – 54.1	38.6	106.9
FAPAS Allergens in Cake Mix - 27364A - July – September 2023	17.6	8.81 – 26.4	14.1	80.1

FAPAS Allergens in Cake Mix - 27364B - July - September 2023	33.6	16.8 - 50.4	28.7	85.4
FAPAS Allergens in Cake Mix - 27331A - July - September 2022	5.66	2.83 - 8.48	5.40	95.4
FAPAS Allergens in Cake Mix - 27331B - July - September 2022	7.16	3.58 - 10.7	7.41	103.5
FAPAS Allergens in Cooked Biscuit - 27314 - January - March 2022	21.8	10.9 - 32.6	18.3	83.9
FAPAS Allergens in Cake Mix - 27311A - November 2021 - February 2022	DETECTED	-	DETECTED	-
FAPAS Allergens in Cake Mix - 27311B - November 2021 - February 2022	DETECTED	-	DETECTED	-
FAPAS Allergens in Cooked Biscuit - 27286 - January - March 2021	13.0	6.5 - 19.5	15.0	115.4
FAPAS Egg in Cake Mix - 27296A - April - June 2021	NOT DETECTED	-	<0.2	-
FAPAS Egg in Cake Mix - 27296B - April - June 2021	15.38	6.43 - 19.3	12.9	119.2
FAPAS Allergens in Cake Mix - 27301A - July - September 2021	9.03	4.51 - 13.50	5.34	59.1
FAPAS Allergens in Cake Mix - 27301B - July - September 2021	6.72	3.36 - 10.10	3.55	52.9

<b>Proficiency Test</b>	<b>Assignment</b>	<b>Result (mg/kg)</b>	<b>Satisfactory results / results agreeing with assignment (%)</b>
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 17 - 1077 - June 2024	DETECTED	DETECTED	100.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 17 - 1177 - June 2024	NOT DETECTED	NOT DETECTED	100.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 17 - 1277 - June 2024	DETECTED	DETECTED	100.0
FAPAS Gluten, Egg and Milk on Swabs (Cotton) 27366E - September - November 2023	DETECTED	DETECTED	97.0
FAPAS Gluten, Egg and Milk on Swabs (Cotton) 27366F - September - November 2023	NOT DETECTED	NOT DETECTED	97.0
FAPAS Egg, Gluten and Milk on Swabs (Cotton) 27333A - September - November 2022	NOT DETECTED	NOT DETECTED	93.0
FAPAS Egg, Gluten and Milk on Swabs (Cotton) 27333B - September - November 2022	DETECTED	DETECTED	97.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 13 - 1077 - June 2022	DETECTED	DETECTED	100.0
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 13 - 1177 - June 2022	DETECTED	DETECTED	100.0



BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 13 - 1277 - June 2022	NOT DETECTED	<0.4	100.0
--	--------------	------	-------