



Proficiency Test	Assigned value (mg/kg)	Result (mg/kg)	Z-score
FAPAS Milk in Infant Breakfast Cereal - 27398A - June - August 2024	NOT DETECTED	<2.5	-
FAPAS Milk in Infant Breakfast Cereal - 27398B - June - August 2024	33.8	40.0	0.70
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula - 19 - 0777 - April 2024	36.2	38.9	0.30
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula - 19 - 0877 - April 2024	55.4	54.6	-0.06
BIPEA Food Allergens - Milk 77C - 18 - Infant Formula - 18 - 0977 - April 2024	NOT DETECTED	<2.5	-



BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 16 - 1077 - January 2024	47.5	31.6	-0.84
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 16 - 1177 - January 2024	NOT DETECTED	<1.0	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 16 - 1277 - January 2024	50.2	35.4	-0.74
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula - 18 - 0777 - November 2023	33.4	29.9	-0.46
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula - 18 - 0877 - November 2023	8.7	8.2	-0.22
BIPEA Food Allergens - Milk 77C - 17 - 0280 Infant Formula - 17 - 0977 - November 2023	NOT DETECTED	<1.0	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 15 - 1077 - June 2023	54.4	41.1	-1.02
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 15 - 1177 - June 2023	NOT DETECTED	<1.0	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 15 - 1277 - June 2023	43.7	32.2	-0.93
BIPEA Allergens in Industrial Process Water 77D - 14 - 0381 Industrial Process Water - 14 - 1077 - January 2023	NOT DETECTED	<1.0	-
BIPEA Allergens in Industrial Process Water 77D - 14 - 0381 Industrial Process Water - 14 - 1177 - January 2023	26.5	32.6	0.55
BIPEA Allergens in Industrial Process Water 77D - 14 - 0381 Industrial Process Water - 14-1277 - January 2023	29.6	33.6	0.39
BIPEA Food Allergens - Milk 77C - Infant Formula - 16 - 0777 - November 2022	15.7	10.7	-1.35
BIPEA Food Allergens - Milk 77C - Infant Formula - 17 - 0877 - November 2022	96.1	47.9	-1.83
BIPEA Food Allergens - Milk 77C - Infant Formula - 15 - 0977 - November 2022	NOT DETECTED	<1.0	-

BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 13 - 1077 - June 2022	NOT DETECTED	<1.0	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 13 - 1177 - June 2022	DETECTED	DETECTED	-
BIPEA Allergens in Industrial Process Water 77D - Industrial Process Water - 13 - 1277 - June 2022	18.8	25.5	0.96

Proficiency Test	Assigned value (mg/kg) according to R-Biopharm Ridascreen Fast Milk R4652	Range (mg/kg) for $z \leq 2$ according to R-Biopharm Ridascreen Fast Milk R4652	Result (mg/kg)	Recovery (%)
FAPAS Milk and Gluten in Infant Soya Formula 27386A February - April 2024	NOT DETECTED	-	<2.5	-
FAPAS Milk and Gluten in Infant Soya Formula 27386B February - April 2024	20.3	10.2 - 30.5	28.2	138.9
FAPAS Allergens in Cooked Biscuit 27382 January - March 2024	25.5	12.8 - 38.3	18.6	72.9
FAPAS Allergens in Cake Mix - 27364A - July - September 2023	36.9	18.5 - 55.4	41.0	111.1
FAPAS Allergens in Cake Mix - 27364B - July - September 2023	66.1	33.0 - 99.1	83.0	125.6
FAPAS Milk in Infant Breakfast Cereal - 27362A - July - August 2023	NOT DETECTED	-	<2.5	-
FAPAS Milk in Infant Breakfast Cereal - 27362B - July - August 2023	29.7	14.8 - 44.5	28.3	95.3
FAPAS Allergens in Cake Mix - 27331A - July - September 2022	22.9	11.5 - 34.4	23.69	103.4

FAPAS Allergens in Cake Mix - 27331B - July - September 2022	11.6	5.80 - 17.4	10.84	93.4
FAPAS Milk in Infant Breakfast Cereal - 27329A - June - August 2022	33.1	16.5 - 49.6	25.47	76.9
FAPAS Milk in Infant Breakfast Cereal - 27329B - June - August 2022	NOT DETECTED	-	<2.5	-
FAPAS Gluten & Milk in Infant Soya Formula - 27316A - February - April 2022	DETECTED	-	DETECTED	-
FAPAS Gluten & Milk in Infant Soya Formula - 27316B - February - April 2022	NOT DETECTED	-	<2.5	-
FAPAS Allergens in Cooked Biscuit - 27314 - January - March 2022	27.7	13.9 - 41.6	31.9	115.2
FAPAS Milk in Infant Soya Formula - 27313A - January - February 2022	NOT DETECTED	-	<2.5	-
FAPAS Milk in Infant Soya Formula - 27313B - January - February 2022	DETECTED	-	DETECTED	-
FAPAS Milk in Infant Breakfast Cereal - 27322A - September - October 2021	NOT DETECTED	-	<2.5	-
FAPAS Milk in Infant Breakfast Cereal - 27322B - September - October 2021	24.3	12.2 - 36.5	24.34	101.4
FAPAS Milk in Infant Formula (hydrolyzed) - 27304A - August - October 2021	NOT DETECTED	-	<2.5	-
FAPAS Milk in Infant Formula (hydrolyzed) - 27304B - August - October 2021	DETECTED	-	DETECTED	-
FAPAS Allergens in Cake Mix - 27301A - July - September 2021	19.6	9.8 - 29.4	14.79	75.5



FAPAS Allergens in Cake Mix - 27301B - July - September 2021	9.78	4.89 - 14.70	7.84	80.2
FAPAS Allergens in Cooked Biscuit - 27286 - January - March 2021	63.6	31.8 - 95.5	48.81	76.8