



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test - 04519 – Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July – September 2024	9.32	7.63	-0.8
FAPAS Food Chemistry Proficiency Test - 04520 - Multi - Mycotoxins in Maize Flour - July – September 2024	6.15	6.69	0.4
FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April – May 2024	6.03	5.82	-0.2



BIPEA 31B - Cereals – Bran - 4 - 3631 – April 2024	4.04	4.68	0.53
BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024	5.83	4.68	-0.67
BIPEA 31B - Mycotoxins – Cereals – Ground Corn - 6 - 187 – December 2023	5.87	8.78	1.69
FAPAS Food Chemistry Proficiency Test - 04497 - Aflatoxins in Peanut Butter (Water Slurry) - November 2023 – January 2024	3.58	3.89	0.4
BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Peanut Paste - 6 - 185 – September 2023	8.69	9.57	0.36
FAPAS Food Chemistry Proficiency Test - 04482 - Multi-Mycotoxins in Maize Flour - July - September 2023	4.72	4.73	0.0
FAPAS Food Chemistry Proficiency Test - 04472 – Multi - Mycotoxins in Maize Flour - April - June 2023	4.73	4.39	-0.3
FAPAS Food Chemistry Proficiency Test - 22199 – Multi - Mycotoxins in Wheat Flour - November 2022- January 2023	3.78	3.34	-0.5
FAPAS Food Chemistry Proficiency Test - 04449 - Mycotoxins in Dried	3.32	3.52	0.3



Figs (Water/Fruit Slurry) - August - September 2022			
FAPAS Food Chemistry Proficiency Test - 04448 – Multi - Mycotoxins in Maize Flour - July - September 2022	3.32	3.71	0.5
FAPAS Food Chemistry Proficiency Test – 04439 – Multi - Mycotoxins in Maize Flour - April - May 2022	5.55	5.60	0.0
FAPAS Food Chemistry Proficiency Test - 04438 - Aflatoxins in Animal Feed (Cereal Based) - April - May 2022	7.80	7.75	0.0
A.I.A. Ring Test - Lotto RTB1 220322 - Aflatoxin B1 in Corn - March 2022	Negative Sample	<limit of detection	-
	9.22	10.17	0.32
	17.76	19.26	0.31
			0.52
FAPAS Food Chemistry Proficiency Test - 22187 – Multi - Mycotoxins in Wheat Flour - November 2021 – January 2022	2.60	2.88	0.5
FAPAS Food Chemistry Proficiency Test - 04426 – Multi - Mycotoxins in Maize Flour - October – December 2021	5.82	6.37	0.3
FAPAS Food Chemistry Proficiency Test - 04418 – Multi - Mycotoxins in Maize Flour - July – September 2021	3.12	3.35	0.3



FAPAS Food Chemistry Proficiency Test - 04417 - Mycotoxins in Hazelnut (Water/Nut Slurry) - June - August 2021	4.20	4.81	0.7
FAPAS Food Chemistry Proficiency Test 04411 Multi-Mycotoxins in Maize Flour April - June 2021	5.35	5.25	-0.1
FAPAS Food Chemistry Proficiency Test - 04408 - Aflatoxins in Peanut (Water/Nut Slurry) - March - April 2021	4.86	5.64	0.7
FAPAS Food Chemistry Proficiency Test - 04402 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2020 - January 2021	20.4	18.18	-0.5
FAPAS Food Chemistry Proficiency Test - 04390 - Mycotoxins in Hazelnut (Water/Nut Slurry) - July - August 2020	4.50	3.67	-0.8
FAPAS Food Chemistry Proficiency Test - 04384 - Multi - Mycotoxins in Maize Flour - April - May 2020	6.13	5.95	-0.1
FAPAS Food Chemistry Proficiency Test - 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	6.65	6.48	-0.1
FAPAS Food Chemistry Proficiency Test - 04375 - Mycotoxin	14.6	12.78	-0.6



Contamination in Animal Feed (Cereal Based) - December 2019 - January 2020			
A.I.A. Ring Test - Lotto RTB1 240919 - Aflatoxin B1 in Corn - September 2019	<limit of detection	<limit of detection	- 0.142
	7.2	7.15	
	13.9	12.90	
	22.3	22.07	