




# BIO-SHIELD FUMONISIN 5

Cat.Number B5548/B5596

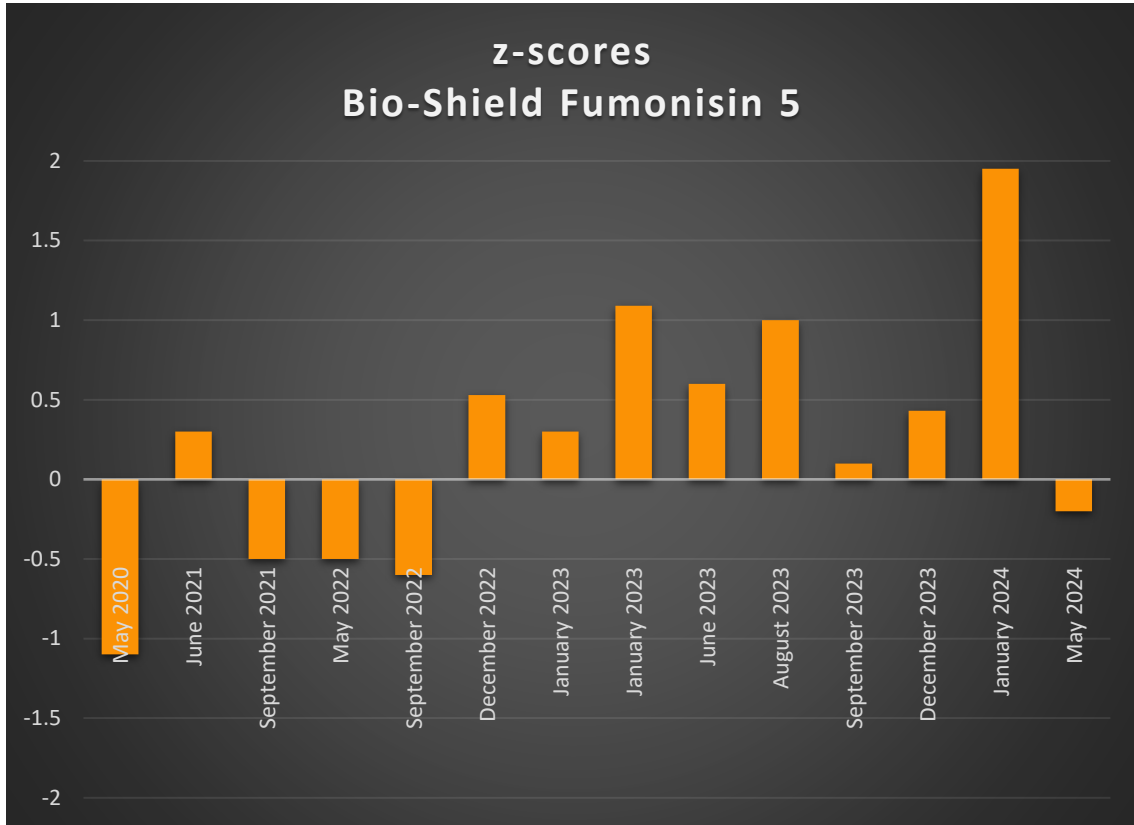

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April - May 2024	788	759.07	-0.2
BIPEA 31E - Mycotoxins - Feed - Animal Feed - 14 - 5 - January 2024	150	226	1.95
BIPEA 31B - Mycotoxins - Cereals - Ground Corn - 6 - 187 - December 2023	434	481	0.43
FAPAS Food Chemistry Proficiency Test - 04482 - Multi - Mycotoxins in Maize Flour - July - September 2023	640	646.3	0.1



## BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04481 - Mycotoxins in Pet Dog Food - July - August 2023	452	533	1.0
FAPAS Food Chemistry Proficiency Test - 04472 - Multi - Mycotoxins in Maize Flour - April - June 2023	798	882	0.6
BIPEA 31E - 3 - Mycotoxin Contamination in Animal feed - 13 - 3031 - January 2023	134	172	1.09
FAPAS Food Chemistry Proficiency Test - 04461 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2022 - January 2023	347	368.87	0.3
BIPEA 31B - 182 - Mycotoxin Contamination in Ground corn - 18 - 0731 - December 2022	167	190	0.53
FAPAS Food Chemistry Proficiency Test - 04448 - Multi - Mycotoxins in Maize Flour - July - September 2022	569	512.5	-0.6
FAPAS Food Chemistry Proficiency Test - 04439 - Multi - Mycotoxins in Maize Flour - April - May 2022	983	911.5	-0.5
FAPAS Food Chemistry Proficiency Test - 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021	720	658.5	-0.5
FAPAS Food Chemistry Proficiency Test - 04411 - Multi - Mycotoxins in Maize Flour - April - June 2021	962	1015.13	0.3



# BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04384 - Multi - Mycotoxins in Maize Flour - April - May 2020	752	612.19	-1.1
--	-----	--------	------

