




BIO-SHIELD OCHRATOXIN 8

Cat.Number B5448/B5496

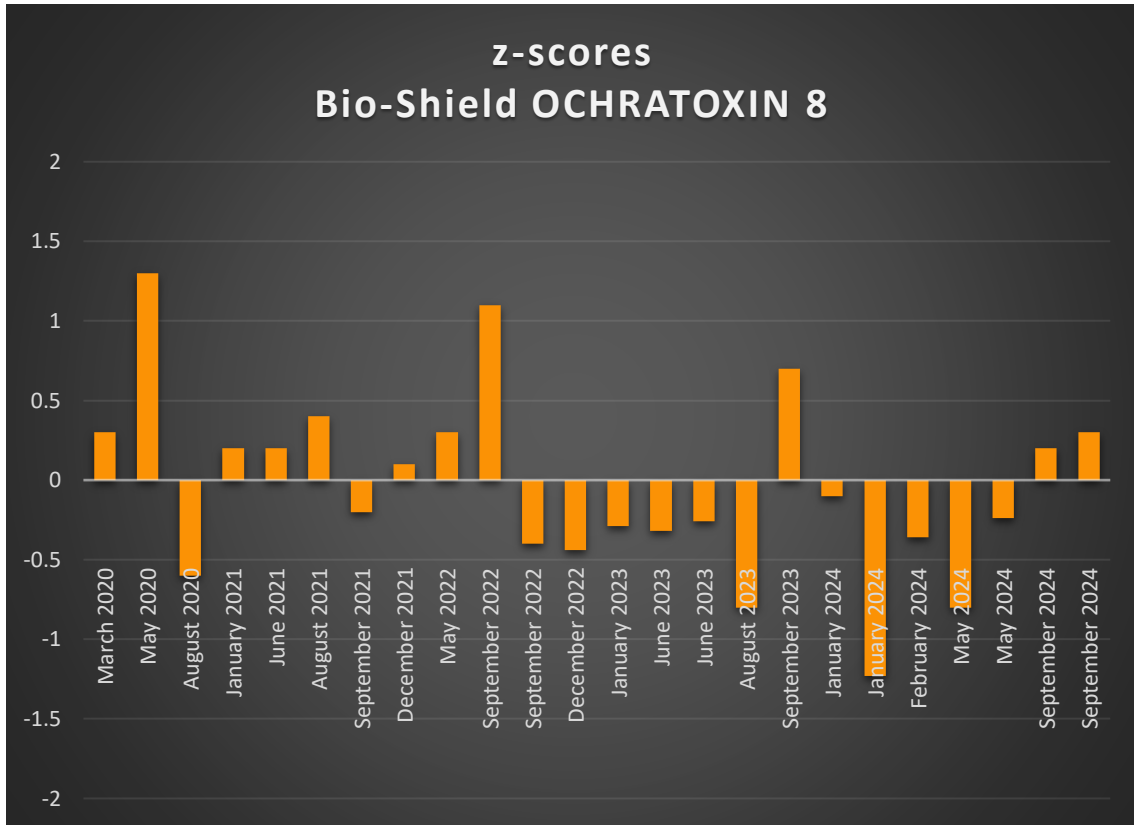

Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
FAPAS Food Chemistry Proficiency Test - 04519 - Multi - Mycotoxins in Pet Dog Food (Dry, Cereal Based) - July - September 2024	15.3	16.4	0.3
FAPAS Food Chemistry Proficiency Test - 04520 - Multi - Mycotoxins in Maize Flour - July - September 2024	2.38	2.47	0.2
BIPEA 31A - Mycotoxins - Dried Fruits, Spices and Other Products - Lentils - 2 - 5931 - May 2024	11.20	10.46	-0.24



BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April - May 2024	4.64	3.82	-0.8
BIPEA 31B - 7 - 188 - Mycotoxins - Cereals - Oats - 7 - 2931 - February 2024	6.65	5.96	-0.36
BIPEA 31E - Mycotoxins - Feed - Animal Feed - 14 - 5 - January 2024	9.31	6.14	-1.23
FAPAS Food Chemistry Proficiency Test - 04498 - Mycotoxin Contamination in Animal Feed (Cereal Based) - November 2023 - January 2024	88.8	86.5	-0.1
BIPEA 31B - Mycotoxins - Cereals - Ground Corn - 6 - 187 - December 2023	Not Set	<1.5	-
FAPAS Food Chemistry Proficiency Test - 04482 - Multi - Mycotoxins in Maize Flour July - September 2023	2.52	2.88	0.7
FAPAS Food Chemistry Proficiency Test - 04481 - Mycotoxins in Pet Dog Food - July - August 2023	10.9	9.13	-0.8
BIPEA 31C - 20 - Mycotoxins - Baby Food - 21 - 3931 - June 2023	2.58	2.38	-0.26
BIPEA 31B - 185 Mycotoxin Contamination in Cereals - Rice - 06 - 4331 - June 2023	4.00	3.62	-0.32
BIPEA 31E - 3 - Mycotoxin Contamination in Animal feed - 13 - 3031 - January 2023	6.07	5.55	-0.29



BIPEA 31B - 183 Mycotoxin Contamination in Ground barley - 05 - 0531 - December 2022	9.08	7.97	-0.44
FAPAS Food Chemistry Proficiency Test - 17227 - Mycotoxins Ochratoxin A in Wine - August - September 2022	2.78	2.40	-0.6
FAPAS Food Chemistry Proficiency Test - 04449 - Mycotoxins in Dried Figs (Water/Fruit Slurry) - August - September 2022	10.4	9.45	-0.4
FAPAS Food Chemistry Proficiency Test - 04448 - Multi - Mycotoxins in Maize Flour - July - September 2022	2.72	3.38	1.1
FAPAS Food Chemistry Proficiency Test - 04439 - Multi - Mycotoxins in Maize Flour - April - May 2022	2.98	3.18	0.3
FAPAS Food Chemistry Proficiency Test - 04426 - Multi - Mycotoxins in Maize Flour - October - December 2021	3.31	3.40	0.1
FAPAS Food Chemistry Proficiency Test - 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021	3.79	3.60	-0.2
FAPAS Food Chemistry Proficiency Test - 04417 - Mycotoxins in Hazelnut (Water/Nut Slurry) - June - August 2021	7.35	7.93	0.4
FAPAS Food Chemistry Proficiency Test - 04411 - Multi - Mycotoxins in Maize Flour - April - June 2021	4.45	4.68	0.2



BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04402 - Mycotoxin Contamination in Animal Feed (Cereal Based) - December 2020 - January 2021	45.7	47.3	0.2
FAPAS Food Chemistry Proficiency Test - 04390 - Mycotoxins in Hazelnut (Water/Nut Slurry) - July - August 2020	7.96	6.87	-0.6
FAPAS Food Chemistry Proficiency Test - 04384 - Multi - Mycotoxins in Maize Flour - April - May 2020	4.82	6.19	1.3
FAPAS Food Chemistry Proficiency Test - 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	6.64	7.06	0.3