



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31B - Mycotoxins - Cereals - Triticale - 3 - 1931 - June 2024	79	70	-0.42
FAPAS Food Chemistry Proficiency Test - 04510 - Multi - Mycotoxins in Maize Flour - April - May 2024	111	97.23	-0.6
BIPEA 31B - Cereals - Bran - 4 - 3631 - April 2024	102	140	1.41
BIPEA 31B - 7 - 188 - Mycotoxins - Cereals - Oats - 7 - 2931 - February 2024	223	209	-0.24



## BIO-SHIELD

BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024	94	85	-0.36
FAPAS Food Chemistry Proficiency Test - 04500 - Multi - Mycotoxins in Durum Wheat - December 2023 – January 2024	80.9	94.0	0.7
BIPEA 31B - Mycotoxins – Cereals – Ground Corn - 6 - 187 – December 2023	210	244	0.63
BIPEA 31B – 186 – Mycotoxin Contamination in Cereals – Common Wheat – 3 - 0631 - October 2023	86.0	64.0	-0.96
FAPAS Food Chemistry Proficiency Test 04482 - Multi-Mycotoxins in Maize Flour - July - September 2023	104	103.2	0.0
FAPAS Food Chemistry Proficiency Test 04481 - Mycotoxins in Pet Dog Food - July – August 2023	40.2	39.1	-0.1
BIPEA 31c 20 Mycotoxins - Baby Food - 21 - 3931 – June 2023	125	129	0.12
BIPEA 31B – 185 Mycotoxin Contamination in Cereals – Rice - 06-4331 - June 2023	139	103	-0.99
FAPAS Food Chemistry Proficiency Test 04472 - Multi-Mycotoxins in Maize Flour - April – June 2023	169	166.7	-0.1
BIPEA 31B-183 Mycotoxin Contamination in Ground barley - 05-0531 - December 2022	84.0	83.0	-0.04



BIPEA 31E-3 Mycotoxin Contamination in Animal feed - 13-3031 - January 2023	113	137	0.8
FAPAS Food Chemistry Proficiency Test 22199 - Multi-Mycotoxins in Wheat Flour - November 2022-January 2023	93.7	72.54	-1.0
BIPEA 31B-182 Mycotoxin Contamination in Ground corn - 18-0731 - December 2022	116	103	-0.43
FAPAS Food Chemistry Proficiency Test 04448 - Multi-Mycotoxins in Maize Flour - July - September 2022	92.6	123.42	1.5
BIPEA 31B-180 Mycotoxin Contamination in Rice - 05-4331 - June 2022	99	66	-1.25
FAPAS Food Chemistry Proficiency Test 04439 - Multi-Mycotoxins in Maize Flour - April - May 2022	189	190.85	0.0
FAPAS Food Chemistry Proficiency Test 22187 - Multi-Mycotoxins in Wheat Flour - November 2021 - January 2022	54	61.1	0.6
FAPAS Food Chemistry Proficiency Test 04418 - Multi - Mycotoxins in Maize Flour - July - September 2021	111	122.64	0.5
FAPAS Food Chemistry Proficiency Test 04411 - Multi-Mycotoxins in Maize Flour - April - June 2021	174	169.47	-0.1
FAPAS Food Chemistry Proficiency Test 22177 - Multi - mycotoxins in Wheat Flour - November 2020 - January 2021	53.5	65.24	1.0



## BIO-SHIELD

FAPAS Food Chemistry Proficiency Test 22175 - Deoxynivalenol, Zearalenone, T-2 and HT-2 in Animal Feed - September - October 2020	523	511.96	-0.1
FAPAS Food Chemistry Proficiency Test 04384 - Multi Mycotoxins in Maize Flour - April - May 2020	330	247.37	-1.3