



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31E - Mycotoxins - Feed - Soyabean Meal - 16 - 6031 - May 2024	12.75	18.26	1.60
BIPEA 31A - Mycotoxins - Dried Fruits, Spices and Other Products - Lentils - 2 - 5931 - May 2024	14.41	12.49	-0.50
BIPEA 31B - Cereals - Bran - 4 - 3631 - April 2024	15.76	19.50	0.89
BIPEA 31B - 7 - 188 - Mycotoxins - Cereals - Oats - 7 - 2931 - February 2024	7.02	6.55	-0.23



## BIO-SHIELD

BIPEA 31E - Mycotoxins – Feed – Animal Feed - 14 - 5 – January 2024	17.86	14.8	-0.65
FAPAS Food Chemistry Proficiency Test - 22215 - Multi - Mycotoxins in Wheat Flour - November 2023 – January 2024	5.90	5.44	-0.4
BIPEA 31B - Mycotoxins – Cereals – Ground Corn - 6 - 187 – December 2023	19.43	15.44	-0.78
FAPAS Food Chemistry Proficiency Test - 04497 - Aflatoxins in Peanut Butter (Water Slurry) - November 2023 – January 2024	8.91	11.07	1.1
FAPAS Food Chemistry Proficiency Test - 04496 - Aflatoxins in Rice – November 2023 – January 2024	9.77	9.02	-0.4
BIPEA 31B – 186 – Mycotoxin Contamination in Cereals – Common Wheat – 3 - 0631 - October 2023	5.06	5.45	0.26
BIPEA 31B – 185 – Mycotoxin Contamination in Cereals – Rice – 06 - 4331 - June 2023	13.56	11.08	-0.68
BIPEA 31B – 184 – Mycotoxin Contamination in Cereals - 02-5431 - April 2023	8.13	8.89	0.33
FAPAS Food Chemistry Proficiency Test - 22199 – Multi - Mycotoxins in Wheat Flour - November 2022 - January 2023	8.88	9.11	0.1
BIPEA 31B – 183 – Mycotoxin Contamination in Ground Barley – 05 - 0531 - December 2022	19.32	20.50	0.23





## BIO-SHIELD

BIPEA 31B - 182 Mycotoxin Contamination in Ground Corn - 18 - 0731 - December 2022	13.30	13.50	0.06
FAPAS Food Chemistry Proficiency Test - 04449 - Mycotoxins in Dried Figs (Water/Fruit Slurry) - August - September 2022	9.34	10.77	0.7
FAPAS Food Chemistry Proficiency Test - 04438 - Aflatoxins in Animal Feed (Cereal Based) - April - May 2022	19.10	16.78	-0.5
FAPAS Food Chemistry Proficiency Test - 22187 - Multi - Mycotoxins in Wheat Flour - November 2021 - January 2022	6.69	6.79	0.1
FAPAS Food Chemistry Proficiency Test - 04417 - Mycotoxins in Hazelnut (Water/Nut Slurry) - June - August 2021	10.2	10.42	0.1
QUALINOVA ROUND : AL 21008 - Aflatoxins in dried fruit - April 2021	9.64	9.07	-0.27
FAPAS Food Chemistry Proficiency Test - 04408 - Aflatoxins in Peanut (Water/Nut Slurry) - March - April 2021	12.1	10.55	-0.6
FAPAS Food Chemistry Proficiency Test - 04390 - Mycotoxins in Hazelnut (Water/Nut Slurry) - July - August 2020	9.67	8.79	-0.4
FAPAS Food Chemistry Proficiency Test - 04378 - Multi - Mycotoxins in Maize Flour - February - March 2020	6.87	6.99	0.1



# BIO-SHIELD

FAPAS Food Chemistry Proficiency Test - 04374 - Aflatoxins in Rice - November 2019 - January 2020	9.31	8.4	-0.4
---	------	-----	------