



Test	Assigned value (µg/kg)	Result (µg/kg)	Z-score
BIPEA 31C - Mycotoxins – Baby Food – 23 - 3931 – June 2024	39.63	24.80	-1.46
BIPEA 31A - 188 - Mycotoxins – Dried Fruits, Spices and Other Products – Hazelnut Paste – 3 - 2031 – March 2024	19.51	20.60	0.21
BIPEA 31A - 186 Mycotoxins - Dried Fruits, Spices and Other Products – Ground Ginger – 02 - 1231 - November 2023	20.96	21.20	0.04

BIO-SHIELD

BIPEA 31A - Mycotoxins – Dried Fruits, Spices and Other Products – Peanut Paste - 6 - 185 – September 2023			8.69	9.57	0.36
BIPEA 31c 20 Mycotoxins - Baby Food - 21 - 3931 – June 2023			14.69	17.7	0.77
BIPEA 31A - 184 Mycotoxins Dried Fruits, Spices and Other Products – Pistachio Paste - 08-1431 May 2023			21.16	17.10	-0.73
QUALINOVA FOOD Mycotoxins in Feed - AL 23003 - April 2023	Sample 1		<1.0	0.989	-
	Sample 2		6.42	6.31	-0.07
BIPEA 31E-3 Mycotoxin Contamination in Animal feed 13 - 3031 January 2023			11.32	9.57	-0.57
BIPEA 31B-180 Mycotoxin Contamination in Rice 05 - 4331 June 2022			9.62	9.37	-0.09
BIPEA 31E-2 Mycotoxin Contamination in Corn Silage 01 -5831 May 2022			6.48	4.29	-1.17
FAPAS Food Chemistry Proficiency Test 04367 Mycotoxins in Dried Figs August - October 2019			10.8	8.1	-1.1
DRRR RVEP 190663 CORN AFLATOXIN 1	Sample 1		9.0	10.1	0.54
	Sample 2		17.9	16.6	-0.12



BIO-SHIELD

FAPAS Food Chemistry Proficiency Test 04364 Mycotoxins in Paprika June - August 2019	10.2	7.9	-1.0
FAPAS Food Chemistry Proficiency Test 04357 Aflatoxins in Peanut March - April 2019	12.40	13.86	0.5
FAPAS Food Chemistry Proficiency Test 04354 Multi Mycotoxins in Maize February - March 2019	4.95	4.78	-0.2
FAPAS Food Chemistry Proficiency Test 04330 Multi Mycotoxins in Maize February - March 2018	4.03	3.63	-0.4
BIPEA 31b 154 Mycotoxins - Cereals - Bran - 02-3631 - April 2017	28.0	21.5	-0.9