



Test	Assigned value (ng/kg)	Result (ng/kg)	Z-score
BIPEA 31D - Mycotoxins – Dairy Food - Contaminants in Dairy Food – 08 - 0431 May 2024	0.089	0.079	-0.38
A.I.A Ring Test Lotto RT M1 260324 Aflatoxin M1 in Lyophilized Milk March 2024	11.95	11.43	-0.21
	21.49	22.43	0.25
	44.34	45.08	0.12
	58.92	58.27	-0.08
BIPEA 31D - Mycotoxins – Dairy Food - Contamination in Milk Powder 15-0331 November 2023	16.4	22.0	1.14

A.I.A Ring Test Lotto RT M1 260923 Aflatoxin M1 in Lyophilized Milk September 2023	15.49	16.20	0.21
	23.92	24.40	0.09
	41.90	41.50	-0.05
	52.78	49.70	-0.44
A.I.A Ring Test Lotto RT M1 280323 Aflatoxin M1 in Lyophilized Milk March 2023	12.21	10.52	-0.18
	21.50	21.24	-0.07
	40.93	40.92	0.00
	53.85	50.53	-0.40
A.I.A Ring Test Lotto RT M1 270922 Aflatoxin M1 in Lyophilized Milk September 2022	10.24	10.70	0.2
	20.74	21.50	0.2
	40.88	42.20	0.2
	54.95	55.90	0.1
A.I.A Ring Test Lotto RT M1 220322 Aflatoxin M1 in Lyophilized Milk March 2022	13.33	9.10	-1.34
	25.47	21.10	-0.93
	48.35	41.60	-0.75
	63.43	58.00	-0.51
A.I.A Ring Test Lotto RT M1 280921 Aflatoxin M1 in Lyophilized Milk September 2021	8.48	10.90	0.92
	15.03	17.39	0.74
	35.62	34.03	-0.28
	46.80	43.88	-0.44
	26.4	26.73	0.047
	19.5	19.39	

A.I.A Ring Test Lotto RT M1 290920 Aflatoxin M1 in Lyophilized Milk September 2020	44.0	44.72	
	50.0	49.41	
A.I.A Ring Test Lotto RT M1 240320 Aflatoxin M1 in Lyophilized Milk March 2020	19.7	20.74	-0.026
	27.1	27.12	
	46.1	45.79	
	56.9	56.49	
A.I.A. Ring Test Lotto RT M1 240919 Aflatoxin M1 in Lyophilized Milk September 2019	13.00	12.61	-0.48
	31.80	28.12	
	50.1	46.25	
	51.1	45.90	
A.I.A. Ring Test Lotto RT M1 260319 Aflatoxin M1 in Lyophilized Milk March 2019	23.72	21.02	-0.69
	53.35	48.91	
	34.63	30.72	
	70.76	66.03	
A.I.A. Ring Test Lotto RT M1 250918 Aflatoxin M1 in Lyophilized Milk September 2018	10.89	8.60	-0.22
	23.05	22.22	
	49.16	49.72	
	58.88	56.96	
FAPAS Proficiency Test 04346 Aflatoxin M1 in Milk Powder October-November 2018	85.9	95.7	0.5
FAPAS Proficiency Test 04331 Aflatoxin M1 in Milk Powder February-April 2018	70.9	70.2	0.0



A.I.A. Ring Test Lotto RT M1 260917 Aflatoxin M1 in Lyophilized Milk September 2017	50.30	47.52	0.2
	18.02	17.84	
	27.28	26.24	
	10.12	9.49	